



Mercury Analyzer

RA-5A

Simple, Powerful & Multi-Configurable
Reducing Vaporization CVAAS
Mercury Analyzer RA-5A



Simple, yet Powerful !


The RA-5A delivers Ultra-High Sensitivity and Excellent Precision while only requiring 5mL of sample. The RA-5A is also Environmentally-friendly through its direct purge design that reduces reagent consumption and waste production by over 80% versus comparable models without any need for compressed gases.



Best Sensitivity, Accuracy


RA-5A has achieved the highest level of Sensitivity and Accuracy with new innovatory-crafted detector optics.

Sample	N	Average	C.V.
0.05 $\mu\text{g/L}$	5回	0.052 $\mu\text{g/L}$	2.2%
0.10 $\mu\text{g/L}$	5回	0.101 $\mu\text{g/L}$	1.7%

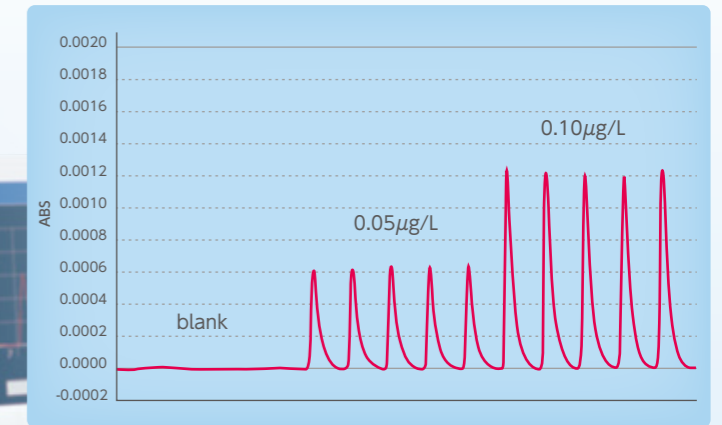

Excellent Repeatability

Memory effect from high mercury concentration is virtually unnoticeable.

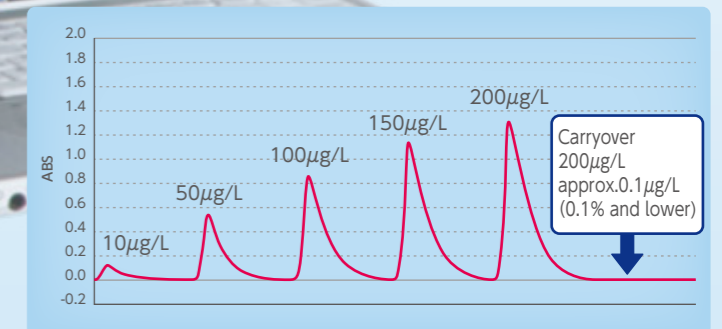

Super Compact

The RA-5A takes up about as much bench space as a sheet of A4-sized paper. Compare to our previous models, both the overall size and weight have been reduced by over 50%. No PC is required for operation, further reducing the occupied bench space.



RA-5A
Setting space is less than 50% (compared to RA-3A)

Options

SC-5 auto sample changer (Sample vessel for 5mL)




SC-5 auto sample changer (Sample vessel for 20mL)



SC-5 auto sample changer (Sample vessel for 100/250mL)



RD-5 auto dispenser

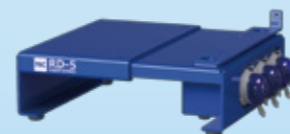


Bottle stand for 5mL (Standard accessory of RD-5)



Bottle stand for 20~250mL



Impinger unit(5mL) (Standard accessory of RA-5A)

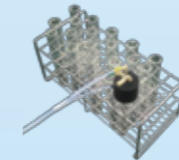


Impinger unit(20mL)



Impinger unit(100/250mL)



SCHEMATIC DIAGRAM

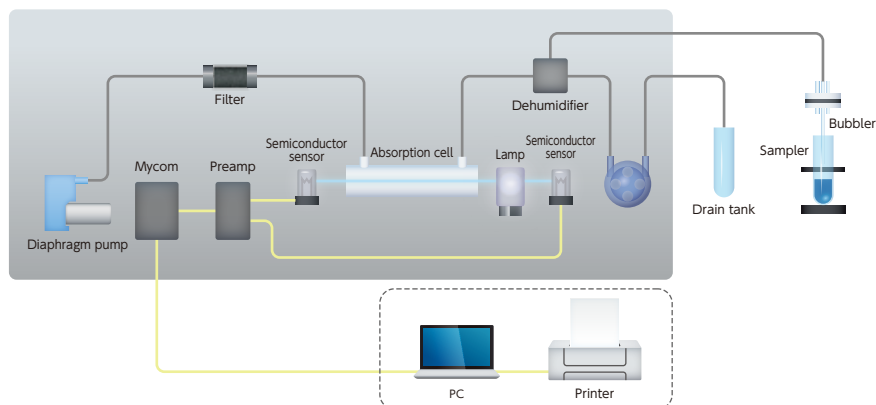

APPLICATIONS

The RA-5a is suitable for a wide variety of aqueous samples, including digested solids. It can be used for mercury analysis in drinking waters, industrial wastewaters, fresh water sources, seawater, soils, sludges, tissues & more.

TEST METHODS

USEPA 245.1, 245.2, 245.5, 7470A, 7471B, ASTM D 3223-02, EN-1483, APHA 3112, JIS K0102 and more

SPECIFICATIONS

Mercury detecting section	Measurement Principle	Non-dispersive double-beam cold vapor atomic absorption spectrometry Patent-pending
	Measurement method	Reducing Vaporization with open air-supply (Peak height measuring and integration measuring selectable)
	Light source	Low pressure mercury discharge lamp
	Wave length	253.7nm
	Detector	Semiconductor sensor (w/254nm bandpass filter)
	Detection Limit	0.0025ng/5mL(0.5ppt)
	Maximum range	2,000ng/5mL(400 ppb)
	Carrier gas required	No Compressed Purified Gas required
	Measuring flow rate	Adjustable (0.1 ~ 0.9 L/min) (just operate with ambient air)
	Measuring time	30sec ~ 999sec
Dehumidifier	Peltier device	
Mercury removal	Removal by activated carbon	
Data processing system (PC-Win / RA-5)	OS	Windows 10 Pro
	Communication	USB
	Display	Peak wave shape, Calibration curve, Measuring time and measuring result
	Data processing software	Selectable from 6 formulas (3 linear and 3 3rd Order fits) , Correlation coefficient is automatically calculated. Statistic calculation : Mean value, standard deviation, Coefficient of variation (All values are automatically calculated.), Self-check function
	Print	Measuring conditions, Calibration curve, Measured values, Statistical calculations, Peak wave shape and Measuring time
	Quality control	Check standard, Avoid abnormally high mercury etc
Installation requirements	Installation dimensions	248W × 253D × 180H (mm)
	Weight	4.7 (kg)
	Installation environment	10°C ~ 35°C, < 10 ~ 80% RH (Non condensing environment is allowed)
Power supply	AC, single phase, 100 ~ 240V 50/60 Hz, 150VA (AC adaptor included)	

SPECIFICATIONS (OPTIONS)

SC-5	Sample vessel : Number	[Standard] : Sample vessel for 5mL : 30 (with vessel sensor function) [Option] : Sample vessel for 20mL : 12 (with vessel sensor function) Sample vessel for 20mL, 100mL and 250mL : 5 (with vessel sensor function)
	Sample access	Random access
	Bubbler tube	PEEK tube (for sample 5mL) / Glass Frit (for sample 20, 100 and 250mL)
	Power supply	AC100 ~ 240V, 50/60Hz, 150VA (AC adaptor included)
	Measurement	180Wx250Dx490H
RD-5	Weight	Approx. 7kg
	Dispensing	Tube pump (3 lines)
	Dispensing volume	Adjustable (Minimum 0 ~ 99.9mL)
	Dispensed reagents	• Sulfuric acid • Tin (II) chloride solution
	Control · Power supply	Direct pump driving from RA-5A
	Measurement	250W x 230D x 80H (with RA-5A connection)
Weight	Approx. 2kg (Stand inclusive)	

STANDARD ACCESSORIES

Sample bottle [Qty 10] / Bottle stand [Qty 1]
/ Impinger [Qty 1] / Bubbler [Qty 2] / Filter unit [Qty 2]

★RA-5A is available to equip with optional attachment accessories, auto sample changer SC-5 and auto dispenser RD-5.
Reagent tube [Qty 1] and reaction tube [Qty 1]
for Optional attachment: SC
Reagent bottle [Qty 1] and Reagent tube [Qty 1]
for Optional attachment :RD

RA-5A series Configurations	RA-5100A	— RA-5A
	RA-5200A	— RA-5A, RD-5
	RA-5300A	— RA-5A, RD-5, SC-5

NIC Nippon Instruments Corporation

Osaka office : 14-8 Akaoji-cho, Takatsuki-shi, Osaka 569-1146 Japan
/Tech. center TEL +81-72-694-5195 FAX +81-72-694-0663
E-mail hg-nic@rigaku.co.jp URL www.hg-nic.com

Singapore office : 10 Science Park Road, #03-24A, Singapore 117684
TEL +65-6873-7068 FAX +65-6873-6372

ISO 9001 : 2015 CERTIFIED Tech-center, Factory & Osaka office/2017. 12. 27
ISO 14001 : 2004 CERTIFIED Tech-center, Factory & Osaka office/2007. 6. 8

Identify the right mercury analyzer for you
<http://www.hg-nic.com>



Distributors

AGS Scientific, Inc.
1 West Bronze Lane
Bryan, TX 77807

Toll Free: 1.877.247.7241
Direct: 979.320.0052
Email: sales@agssci.com
Web: www.agssci.com